

## EzWay™ Coating Buffer (pH 9.6)

**1. Catalog No.** K0331021

**2. Quantity** 125 ml x 2

**3. Storage** Store at 2-8°C. Do not freeze.

**4. Description** ■ Coating buffer for ELISA assay.

**5. Contents**

| Description             | Amount     |
|-------------------------|------------|
| Coating Buffer (pH 9.6) | 125 ml x 2 |

**6. Coating**

- Coating Solution: Resolve the coating material (antigen or antibody) in the coating buffer to make 1 ug/ml (1-10 ug/ml).
- Dispense 100 ul (50-200 ul) of prepared Coating Solution to each well.
- Incubate for overnight at 4°C.