

EzWay™ TMB Solution

- 1. Catalog No.** K0331061
- 2. Quantity** 1000 wells
- 3. Storage** Store at 2-8°C. Do not freeze.

- 4. Description**
- One-step TMB
 - More higher activity level than normal TMB solution

5. Kit Contents

Description	Amount
TMB Solution	100 ml

7. Color Reaction and Reading

- Dispense 100 ul EzWay TMB Solution into each well.
- After sufficient color development (5-30minutes at room temperature or at 37°C), add 100 ul Stop Solution (2M H₂SO₄ or 0.2M H₂SO₄) to each well.
- Read plates in a microwell plate reader at wavelength setting of 450nm